## Dos and Don'ts in the TGADTA and DSC Lab (SAIF IIT Bombay)

- 1. I will abstain from utilizing any personal electronic devices during my designated TA duty hours.
- I will abstain from utilizing any personal returns.
  I will attentively listen to and follow the instructions provided by Meenal and Dr. Nutan while performing my duties.
- 3. I will utilize laboratory instruments with great care and concentration.
- 4. When operating the TGDTA/DSC instrument, I will carefully observe and adhere to the instructions listed on the Material Safety Data Sheet (MSDS), including the heating program and decomposition temperature.
- 5. In DSC, the final heating temperature (as per the heating program) will not surpass the decomposition temperature mentioned on the MSDS.
- 6. Prior to running any samples, I will ensure that the gas supply is properly regulated.
- 7. I will never operate the TGDTA or DSC instrument without turning on the purge gas.
- 8. I will always use appropriate sample pans that are compatible with the instruments.
- 9. Aluminium pans will be used for heating upto a max of 600 °C. Aluminium melts at 660 °C. For heating above 600 °C, I will use either Platinum (Pt) pans or Ceramic pans, and for temperature above 1000 °C, I will use ONLY Ceramic pans.
- 10. I can clearly differentiate between Pt and Aluminium pans. Further I also know that Aluminium pans are one-time use, while Pt pans are very expensive and should be used carefully.
- 11. During DSC runs, I will intermittently monitor the thermogram for any disturbances, such as a continuous wavy baseline, and immediately halt heating or notify Meenal/Dr. Nutan to prevent further sample decomposition, which could damage the DSC cell.
- 12. I will meticulously measure the weights of samples and correctly enter them in the method editor.
- 13. I will handle the weighing balance with extreme care.
- 14. In DSC, the sample pans will be crimped securely to prevent any spills during the runs.
- 15. Plotting of data will be performed with diligence and with a thorough understanding of the specific requirements of the User.
- 16. I will ensure that there is enough Liq. Nitrogen in the Dewar before starting the experiment.
- 17. I will not leave the laboratory before my designated time without notifying.
- 18. Unless due to legitimate health concerns, I will not request leave or any modifications to my duty schedule.
- I will arrive promptly at the designated time for TA duty (9:30 am for forenoon session & 2 pm for the afternoon session).
- 20. I will document all work and samples performed each day in the laboratory log book.
- 21. At the conclusion of each TA duty day, I will record my attendance in the register.